

- Turn the coarse adjustment knob clockwise until the objective is in operating position. You will feel a definite position stop for the objective. Turn the objective clockwise until you are familiar with this position stop.
 - Open the iris diaphragm approximately half way.
 - Place the slide in the stage slide holder securely.
 - View through the eyepieces and without disturbing the coarse-adjustment setting.
- Proper Care of the Microscope:**
- Use both hands when carrying the microscope, one firmly grasping the arm of the microscope, the other beneath the base. Avoid sudden jolts.
 - To keep the microscope and lens systems clean.
 - Never touch the lenses. If the lenses become dirty, wipe them gently with lens paper.
 - Never leave a slide on the microscope when it is not in use.

ADJUST THE BRIGHTNESS

↓

OBSERVE SPECIMEN

↓

RESULT

